

Are you a dairy farmer?

Manage the antibiotic risk
with confidence and at the farm



Detection of beta-lactams (including cefalexin)
and tetracyclins in cow milk

Easy • Fast • Reliable

Adapted
to use
at the farm

WE MANUFACTURE TESTS FOR THE DETECTION OF ANTIBIOTICS AT THE FARM

DipSensor[®]

DETECTION OF BETA-LACTAMS (INCLUDING CEFALEXIN) AND TETRACYCLINES IN COW'S MILK

01



Immerse the tube
in the milk

02




Place the test strip in the tube
in contact with the milk

03



Incubate 10 minutes
at room temperature

VISUAL INTERPRETATION OF THE RESULTS

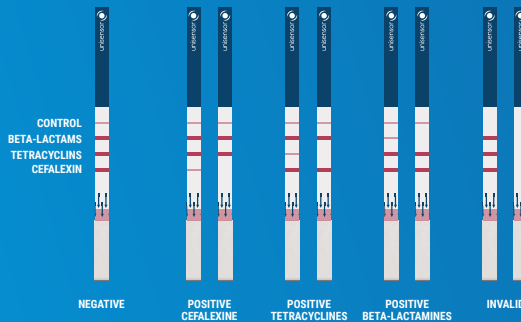

CONTACT US

milktest@intermed.be

International Medical Products
Waversessesteenweg 1110 - 1160 Brussel
Chaussée de Wavre 1110 - 1160 Bruxelles
02/660.50.75 - www.intermed.be