

## Rapid Test Detecting Betalactams, Sulfonamides And Tetracyclins In Milk

TriSensor is a rapid assay in dipstick format detecting the contamination of milk samples by Betalactams, Tetracyclins and Sulfonamides molecules.


The test can be used in-situ for daily controls or in laboratories for the analysis of series of milk samples. This test does not require any sample processing, cleaning or extraction.



### BENEFITS

- X3 MULTIPLE**  
Detection of  $\beta$ -lactams, Tetracyclins and Sulfonamides.
- RAPID**  
Results in 6 minutes.
- SENSITIVE**  
EU MRLS.
- RELIABLE AND ROBUST**  
KG
- USER-FRIENDLY**  
Easily performed on site or in lab.
- COST-EFFECTIVE**  
€

### HOW TO USE

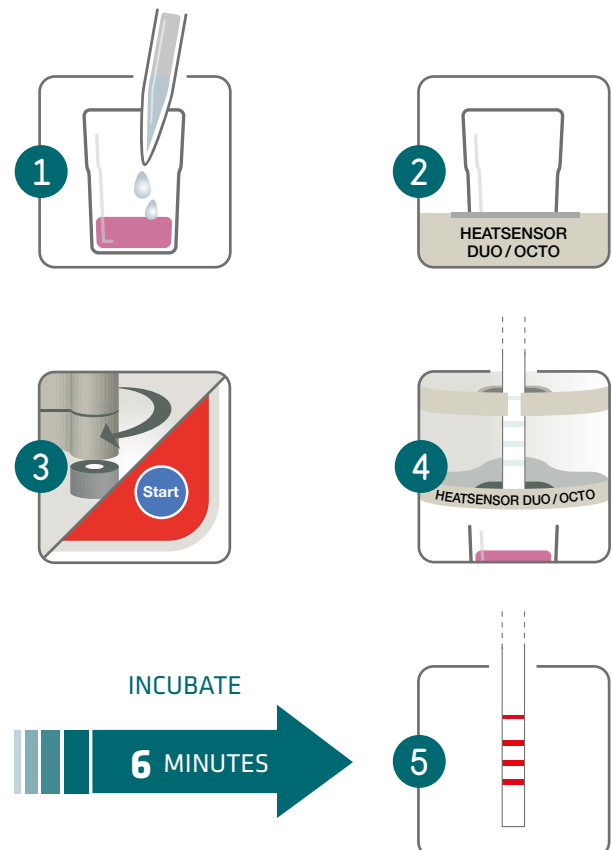


## SENSITIVITY

	<b>β-LACTAMS</b>	<b>LOD (PPB)</b>	<b>SULFONAMIDES</b>	<b>LOD (PPB)</b>	<b>TETRACYCLINES</b>	<b>LOD (PPB)</b>
<b>PENICILLINS</b>	Penicillin-G	2,5-3,5	Sulfadiazine	8-10	Tetracycline (*)	80-100
	Ampicillin	3-4	Sulfapyridine	0,5-1	Oxytetracycline	60-70
	Amoxicillin	3-4	Sulfathiazole	7,5-8,5	Chlortetracycline	50-60
	Oxacillin	13-15	Sulfamethoxazole	320-360	Doxycycline	20-30
	Cloxacillin	7-9	Sulfamethazine	1-2		
	Dicloxacillin	4-6	Sulfamethoxy-pyridazine	2-3		
	Nafcillin	50-70	Sulfadimethoxine	10-15		
<b>CEFALOSPORINS</b>	Ceftiofur	8-12	Sulfacetamide	300-600		
	Cefquinome	16-20	Sulfamerazine	2-3		
	Cefazolin	15-19	Sulfamonomethoxine	8-12		
	Cefapirin	4-6	Sulfaquinoxaline	14-18		
	Cefacetrile	16-20	Sulfachloropyridazine	5-10		
	Cefoperazone	3-4	Sulfaguanidine	15-25		
	Cefalexin	700-800	Sulfamethizole	220-260		
	Cefalonium	3-5	Sulfasalazine	250-350		

(\*) Other Tetracyclines compounds are detected at a lower concentration than Tetracycline: Minocycline, Demeclocycline, Sancycline, Amicycline, Meclocycline.

## RESULTS

Visual interpretation of the result obtained by comparing the intensity of each test line with the control (CTRL) line.

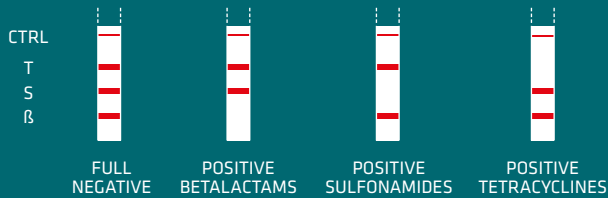

T: Tetracyclines - S: Sulfamides - β: Betalactams

An instrumental reading is also possible with the ReadSensor 2 device.

# TriSensor

## PRODUCTS REFERENCES

**KIT035**  
TRISENSOR MK BTS MRL - 96 PC

**APP032**  
HEATSENSOR DUO

**APP052**  
HEATSENSOR OCTO



**APP088**  
READSENSOR 2



**unisensor**  
DIAGNOSTIC ENGINEERING

Liège Science Park  
Rue Louis Plescia, 8  
B-4102 Ougrée (Seraing)  
BELGIUM

[www.unisensor.be](http://www.unisensor.be)

Phone +32 4 252 66 02  
[info@unisensor.be](mailto:info@unisensor.be)

Reseller

**International Medical Products S.A.**  
Marcel Broodthaersplein 8/5 Place Marcel Broodthaers  
BRUSSEL 1060 BRUXELLES - BELGIUM  
T: +32 (0)2 660 50 75 - F: +32 (0)2 660 20 98  
[www.intermed.be](http://www.intermed.be) - [info@intermed.be](mailto:info@intermed.be)

**beldico** B.V.  
Verlengde Poolseweg 16  
4818 CL Breda  
T: +31 (0) 765 244 663  
[www.beldico.nl](http://www.beldico.nl) - [info@beldico.nl](mailto:info@beldico.nl)